

Wear eye protection.

Wear gloves.

Consult SDS for proper PPE.



1



Using gloves and eye protection, insert a dry roll of Sanotex<sup>®</sup> wipes into the Sanotex<sup>®</sup> canister.

NOTE: For 10" x 12" rolls (#85340 or #85340), place the roll with folded wipe edge up and the unfolded edge down.

2



Carefully pour disinfectant solution over center of wipe roll in canister as noted below. Replace lid.

Wait 5 minutes before use to assure proper saturation of wipes.

Part Number	Description	Amount of Disinfectant Solution
85338	10 x 12" (40 ct.)	16 oz.
85339	6 x 8" (150 ct.)	24 oz.
85340	10 x 12" (60 ct.)	24 oz.

3



After 5 minutes, remove the lid. From the center of the roll, pull a wipe up and carefully thread it through the opening in the Sanotex canister lid approximately one inch.

Replace lid, screwing on tightly.

Apply secondary label and write date on which wipes were saturated with a permanent marker.

4



To dispense, open the hinged lid and pull a single, saturated wipe through the opening at an angle. The spring-loaded lid will re-close automatically.

Always wear gloves and eye protection when preparing and using saturated wipes.

### Questions?

Call Contec at 1-800-289-5762 or your distributor at \_\_\_\_\_